MISTRAL EVOLUTION NITROGEN GENERATOR





North America & South America contact:

Europe, Asia, Africa, & Australia/Oceania contact:

VICI DBS USA

tel: +1 713-263-6970 fax: +1 713-263-6971 web: www.vicidbs.com

VICI AG INTERNATIONAL

tel: +41 41 925-6200 fax: +41 41 925-6201 web: www.vicidbs.com

DESCRIPTION

The VICI DBS° Mistral Evolution is a self contained generator that produces up to 40 L/min of pure LC/MS grade nitrogen at up to 8 barg (116 psig). Nitrogen is produced by utilizing a combination of compressor and Carbon Molecular Sieve (CMS) technology. High and low pressure compressors are carefully matched to the CMS demand to ensure quite and reliable operation. This unique combination of dual compressor technology has several unique advantages over all other nitrogen generators commercially available.

Nitrogen is produced at low pressure, which ensures a longer compressor life and then compressed to 8 barg (116 psig) using a second stage compressor. This combination guarantees a long compressor life reducing maintenance costs and down time.



INCREASE EFFICIENCY

The relatively high gas volumes required by LCMS instruments make cylinder supply inappropriate and liquid nitrogen expensive. A constant, uninterrupted gas supply eliminates interruptions of analyses to change cylinders.



RETURN ON INVESTMENT

Payback period can be as short as 6 to 12 months.



IMPROVE SAFETY

Nitrogen produced at low pressure and ambient temperature removes the hazards associated with high pressure cylinders and liquid dewars.



ENHANCE PERFORMANCE

Production of a constant flow and pressure of nitrogen improves the consistency of the LC/MS analysis, results and reproducibility.



SUPERIOR TECHNOLOGY

The 2-stage pressure design allows the compressors to work at their optimum pressure range reducing the stress and results in extending the life time of the compressors. This reduces maintenance costs and associated downtime.





FEATURES

Produces a continuous supply of high purity nitrogen | On-demand supply 24/7 | Flow rate at 8 barg (116 psig): 8 to 40 L/min | Purity: LC/MS grade | Integrated low noise oil free compressors | Proprietary carbon molecular sieve technology | 2-year complete product warranty | Complete "plug and play" LC/MS solution



BENEFITS

Eliminates dangerous high pressure cylinders helping to keep your employees safer | Removes the logistics, inconvenience, downtime and costs of a cylinder system | Flow capacity to match your specific instrument demands | Ideal for all LC/MS applications | Simple installation and operation | Superior gas purification | Peace of mind | Install directly in the laboratory



APPLICATIONS

LC/MS INSTRUMENTS

- Nebulizing gas
- Curtain gas
- Shield gas
- Sheath gas
- Electrospray gas
- APCI gas
- Jet stream gradient

SPECTROSCOPY

Nuclear magnetic resonance spectrometers (NMR)

ANALYZERS

Thermal analyzers (TGA/DSC)

OTHER APPLICATIONS


- Sample concentrators
- Evaporative light scattering detector (ELSD)

- Corona detectors
- Charged aerosol (CAD)
- Sample evaporators
- Cell incubators
- IVF Stem cell incubators
- Mask aligner
- Fume hood/cabinet purge
- 3D printer inerting
- Glove box purge

OPERATING DIAGRAM

The Mistral Evolution (PSA technology) includes two dynamically balanced, long-life compressors and delivers a continuous or on demand stream of pure nitrogen gas of 99.5% with a flow rate to 40 L/min.

It uses pressure swing adsorption (PSA) system which removes oxygen, carbon dioxide and water from compressed air. The resulting stream of pure nitrogen is ideal for LCMS instruments and other laboratory applications where inert gases are required.

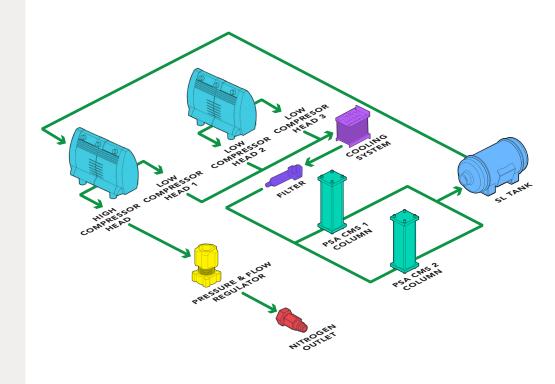

ORDERING INFORMATION (for best service, please call to discuss your application before placing your order).

MISTRAL EVOLUTION 10

DB-EVO-10-EU 220V/50Hz **DB-EVO-10-US** 115V/60Hz

MISTRAL EVOLUTION 25

DB-EVO-25-EU 220V/50Hz **DB-EVO-25-US** 115V/60Hz

MISTRAL EVOLUTION 35

DB-EVO-35-EU 220V/50Hz **DB-EVO-35-US** 115V/60Hz

MISTRAL EVOLUTION 40

DB-EVO-40-EU 220V/50Hz **DB-EVO-40-US** 115V/60Hz